



⚠ Weight Capacity: 50 lbs. / sq. ft. (244kg / m²)

⚠ Worktable is not intended for over-the-road use.

1. Assemble posts.

<p>all mobile units</p> <p>bumper x4</p>	<p>all mobile units with 3-sided bottom frame</p> <p>collar adaptor x4</p> <p>attach to lowest possible location</p>	<p>all stationary units with 3-sided bottom frame</p> <p>x4</p> <p>6" / 152mm</p>
--	--	---

2. Attach 3-sided frame or bottom shelf to posts

Be sure the bumpers are pointing out.
Attach frame or shelf as shown below. Seat the shelf onto the posts with a mallet or a hammer and board.

mobile unit with solid shelf

Note: refer to the manual inside the shelf carton for MetroMax assembly information



mobile unit with 3-sided frame

stationary units
attach to lowest practical location

3. Attach top frame and corner brace supports

Note: corner braces are not included with 30" x 30" (762mm x 762mm)

First, slide the two corner brace supports down over the rear posts. Attach collar adaptors to the tops of all four posts. Attach top frame to posts and collar adaptors.



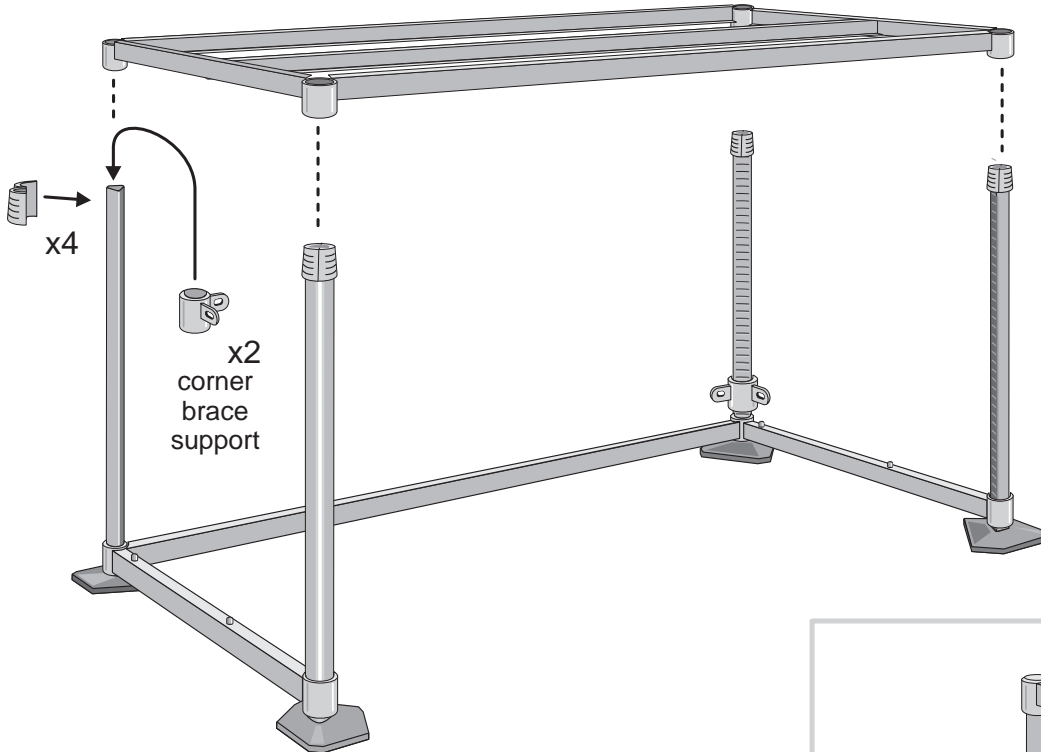
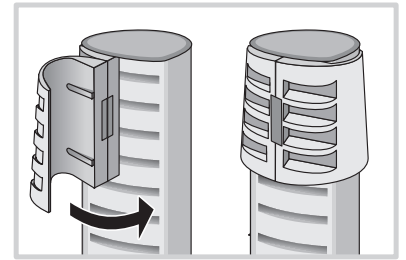


4. Seat top frame

Use a mallet or a hammer and board as shown below.



5. Attach corner braces (except 30" x 30" [762mm x 762mm] unit)

For each of the the two rear posts, lift the corner brace support up and attach a collar adaptor to the post so that when the the support is lowered into place the measurement between the collars is as shown at right. Attach the corner braces as shown using nylok nuts.

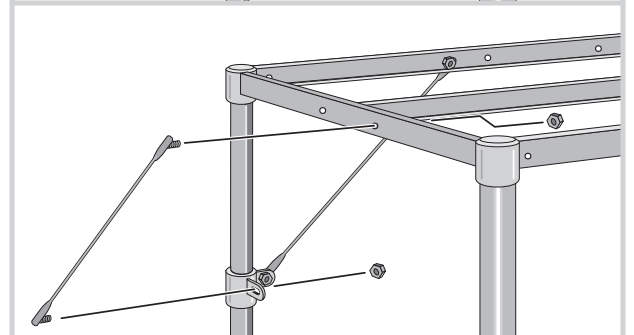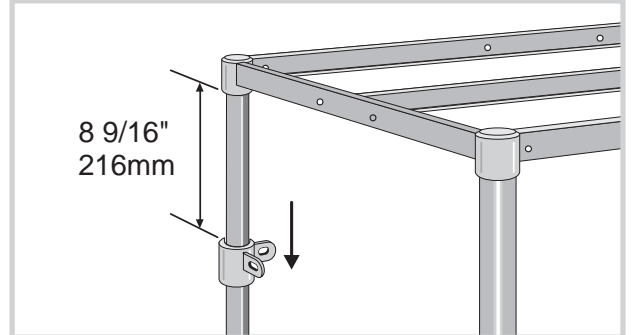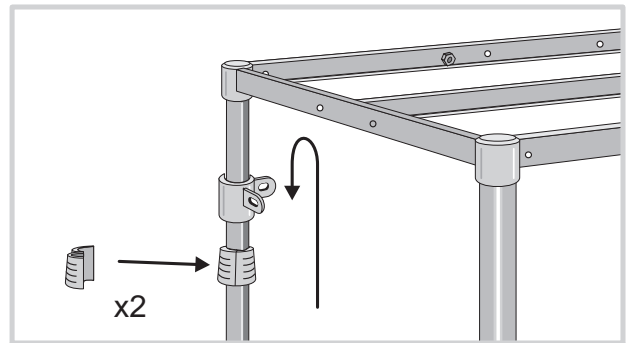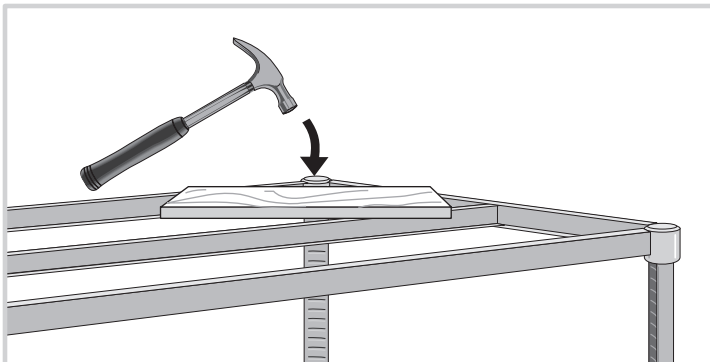


6. Attach optional drawer slides or keyboard tray

Refer to the chart at right for mounting locations. Attach the accessories to the underside of the top frame as shown below. Do not install drawer until the work surface is attached. Lock the keyboard tray in place before attaching the work surface.

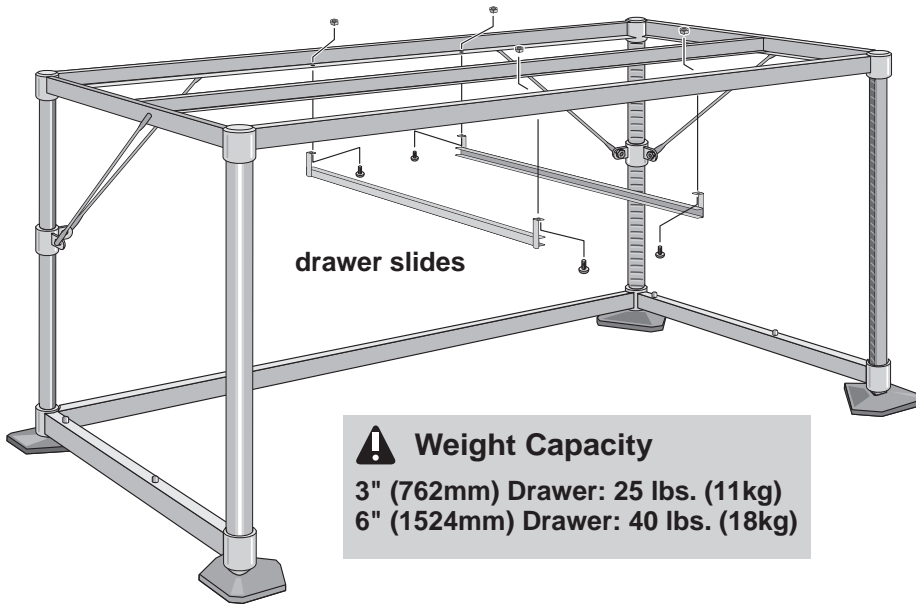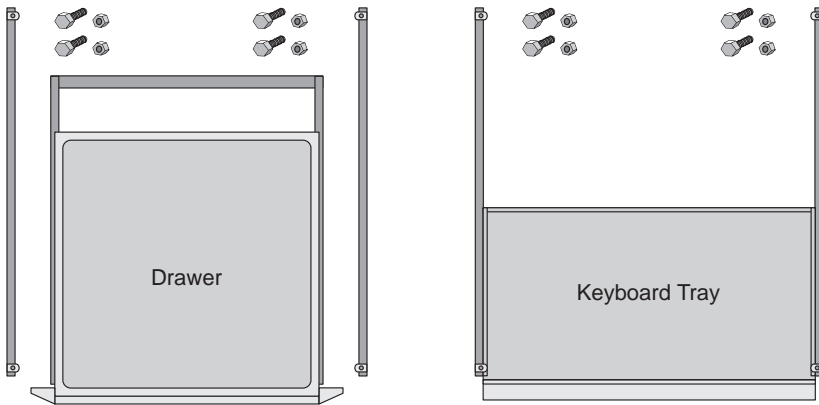

⚠ Weight Capacity
 3" (762mm) Drawer: 25 lbs. (11kg)
 6" (1524mm) Drawer: 40 lbs. (18kg)

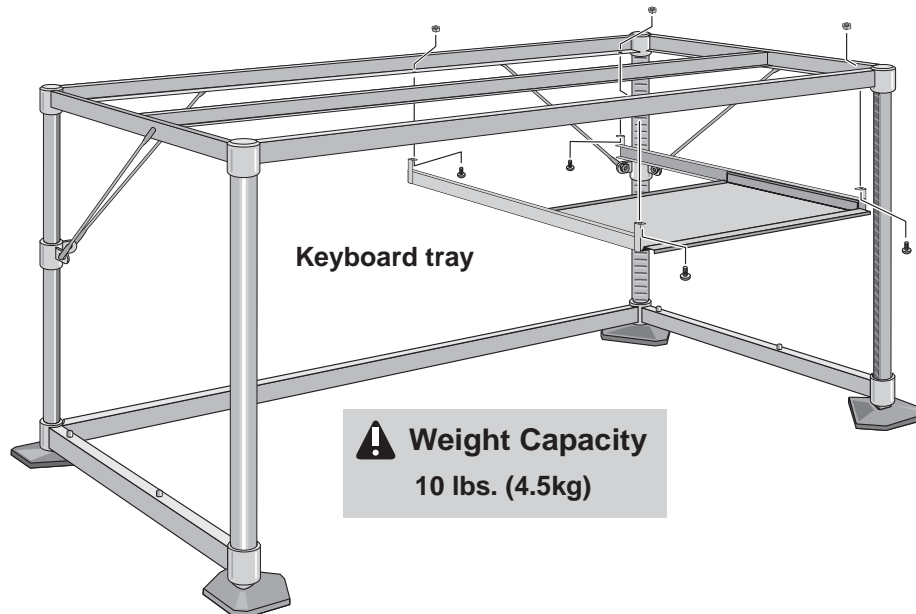

⚠ Weight Capacity
 10 lbs. (4.5kg)

Possible Accessory layouts:

Note: Drawers and the Keyboard Tray may be installed to face either the front or the rear of the table



Drawer




Keyboard Tray

30" x 30"
(760mm x 760mm)


30" x 36"
(760mm x 914mm)



30" x 48" (760mm x 1219mm)

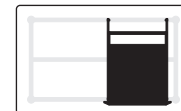

30" x 60" (760mm x 1524mm)

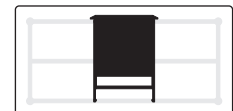


7. Attach worksurface

Place the worksurface upside down on a non-scratching surface on the floor. Identify the rear of the phenolic worksurface by the partially drilled holes along the back edge (for attaching the optional backsplash).

Clean the bottom of the worksurface with household cleaner and remove the backing from the mounting tape on the top frame.

Turn the table frame upside down and carefully lower the frame onto the worksurface. As shown below, align the guidelines on the bottom of the worksurface with the inside edges of the frame. Press the frame down to assure good adhesion.

If the table is mobile, attach the casters (casters were ordered separately).

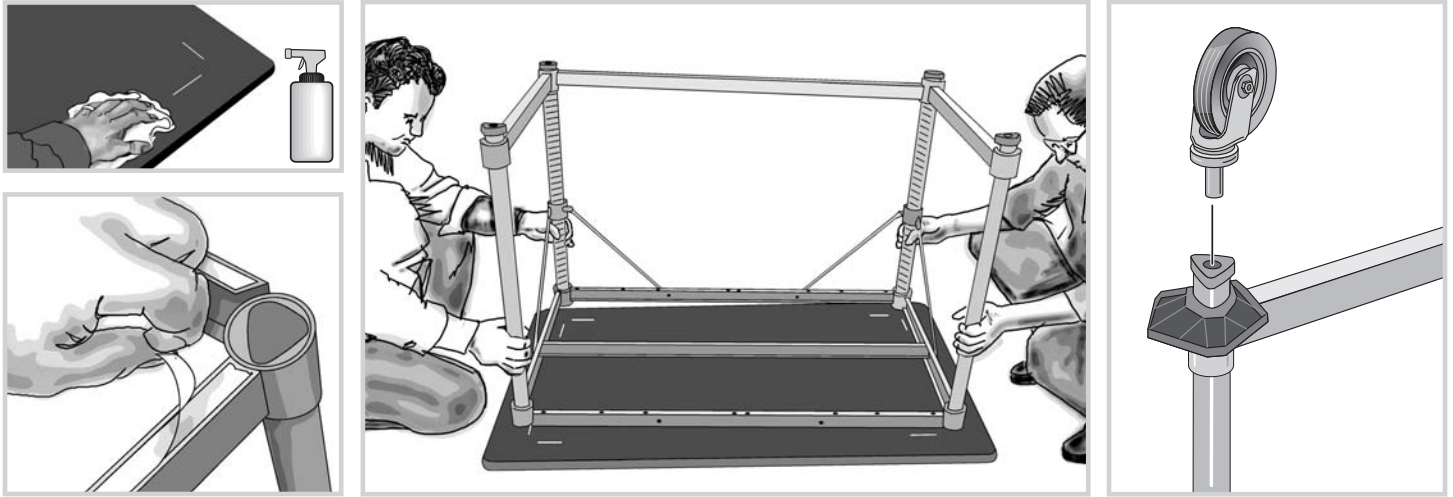


8. Attach optional backsplash (phenolic worksurface)

Remove the protective plastic from the top of the worksurface and the front of the backsplash. Use a drill to drive the drill-point screws up through the partially drilled holes until they break the top surface.

Place the backsplash into position, aligning the holes. Drive the screws up into the holes in the backsplash.

You may optionally apply silicone sealant to the joint between the backsplash and the worksurface.

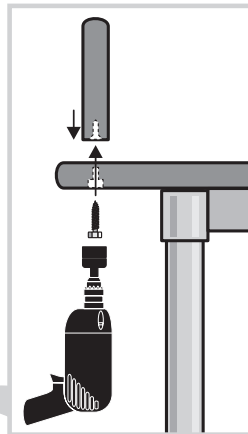

9. Install optional wire shelf onto 3-sided frame

Position the wire shelf over the studs on the 3-sided frame.

⚠ Weight Capacity
150 lbs. (68kg)



InterMetro Industries Corporation

North Washington Street, Wilkes-Barre, PA USA 18705

For Product Information, Call 1-800-433-2232

Pour de plus amples renseignements sur les produits,
appeler au 570-825-2741

Para obtener información sobre el producto llame al: 570-825-2741

Visit our Web Site / Visitez notre site Web / Visite nuestro sitio en la Web: www.metro.com

INSTR-LABTABL
5/05

Information and specifications are subject to change without notice. Please confirm at time of order.
Les données et caractéristiques techniques sont susceptibles d'être modifiées sans préavis. Veuillez confirmer au moment de la commande.
Las informaciones y especificaciones están sujetas a cambios sin previo aviso. Por favor, confirme a la vez que realiza el pedido.